# NVIGATOR™ 系列

## NV/NVT

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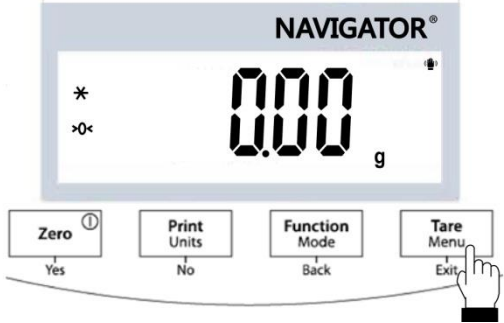
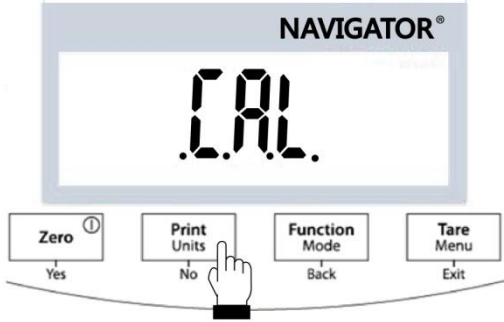
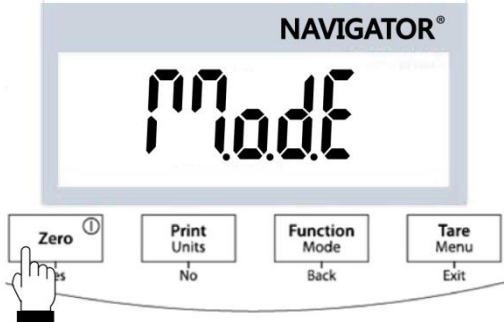
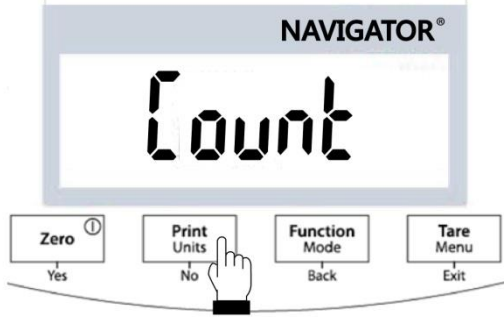
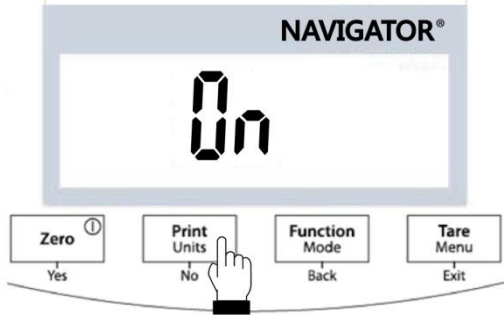
<p>1. 長按 Menu 進入校正鍵，顯示 CAL</p>	 <p>The display shows 'NAVIGATOR' at the top, a '*' symbol on the left, and '0.00 g' in the center. Below the display are four buttons: 'Zero' (with a circled '1'), 'Print Units', 'Function Mode', and 'Tare Menu'. Underneath each button are the options 'Yes', 'No', 'Back', and 'Exit' respectively. A hand icon is shown pressing the 'Tare Menu' button.</p>
<p>2. 顯示 CAL 後按 NO 鍵選擇 Mode 模式</p>	 <p>The display shows 'NAVIGATOR' at the top and 'CAL' in the center. Below the display are the same four buttons as in the previous step. A hand icon is shown pressing the 'Print Units' button.</p>
<p>3. 顯示 Mode 後按 Yes 進入選擇預模式</p>	 <p>The display shows 'NAVIGATOR' at the top and 'Mode' in the center. Below the display are the same four buttons. A hand icon is shown pressing the 'Zero' button.</p>
<p>4. 選擇預開啟或關閉的模式按 Yes 進入</p>	 <p>The display shows 'NAVIGATOR' at the top and 'Count' in the center. Below the display are the same four buttons. A hand icon is shown pressing the 'Print Units' button.</p>
<p>5. 按 NO 選擇開或關，並按 YES 確認</p>	 <p>The display shows 'NAVIGATOR' at the top and 'On' in the center. Below the display are the same four buttons. A hand icon is shown pressing the 'Print Units' button.</p>

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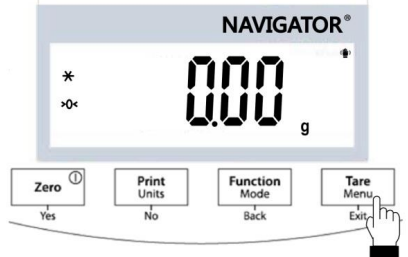
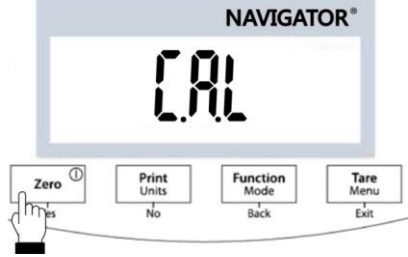
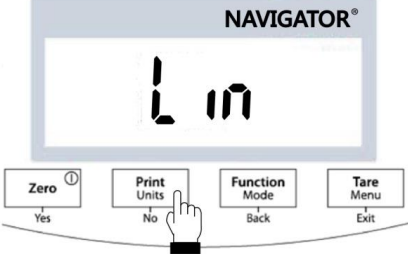
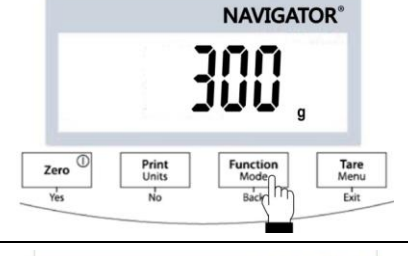
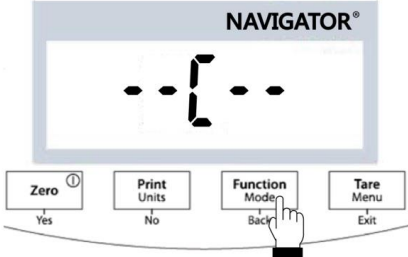
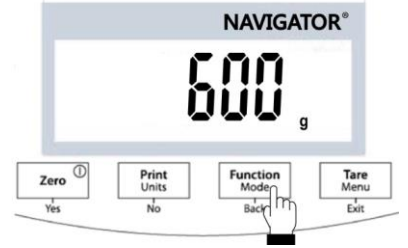
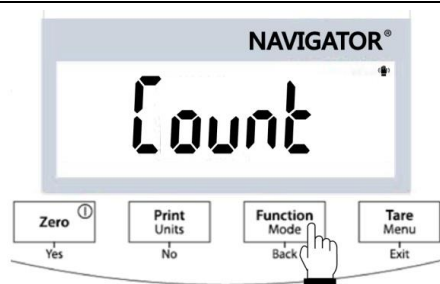
<p>1. 按 <b>Yes</b> 確認，螢幕出現 CLR.RPU</p>	 <p>The screenshot shows the NAVIGATOR scale display with 'CLR.RPU' on the screen. Below the screen are four buttons: 'Zero' (with a circled '1' above it), 'Print Units', 'Function Mode', and 'Tare Menu'. Under 'Zero' are 'Yes' and 'No' options. Under 'Print Units' is 'No'. Under 'Function Mode' is 'Back'. Under 'Tare Menu' is 'Exit'. A hand icon is pointing to the 'Zero' button.</p>
<p>2. 按 <b>Yes</b> 確認，螢幕出現 <b>Pwt 10</b> 訊息而且數字 10。可以按 <b>No</b> 或 <b>Back</b> 來增加或減少數值。 例如，增加數值到 20，請按 <b>No</b>。</p>	 <p>The screenshot shows the NAVIGATOR scale display with 'Pwt 10' and 'PC' on the screen. Below the screen are four buttons: 'Zero' (with a circled '1' above it), 'Print Units', 'Function Mode', and 'Tare Menu'. Under 'Zero' are 'Yes' and 'No' options. Under 'Print Units' is 'No'. Under 'Function Mode' is 'Back'. Under 'Tare Menu' is 'Exit'. A hand icon is pointing to the 'Function Mode' button.</p>
<p>3. 放置 10 個樣品在稱盤上。按 <b>Yes</b> 鍵會以此 10 個樣品來建立平均單重 (APW)。螢幕顯示 10 個(pcs)。</p>	 <p>The screenshot shows the NAVIGATOR scale display with '10' and 'PC' on the screen. Below the screen are four buttons: 'Zero' (with a circled '1' above it), 'Print Units', 'Function Mode', and 'Tare Menu'. Under 'Zero' are 'Yes' and 'No' options. Under 'Print Units' is 'No'. Under 'Function Mode' is 'Back'. Under 'Tare Menu' is 'Exit'. A hand icon is pointing to the 'Zero' button.</p>
<p>4. 從稱盤上移除 10 個樣品，再放置更多的樣品到盤上。螢幕會顯示相對應的數量。</p>	 <p>The screenshot shows the NAVIGATOR scale display with '20' and 'PC' on the screen. Below the screen are four buttons: 'Zero' (with a circled '1' above it), 'Print Units', 'Function Mode', and 'Tare Menu'. Under 'Zero' are 'Yes' and 'No' options. Under 'Print Units' is 'No'. Under 'Function Mode' is 'Back'. Under 'Tare Menu' is 'Exit'. A hand icon is pointing to the 'Zero' button.</p>
<p>5. 要檢視平均單重或樣品數量，按 <b>Function/Mode</b> 按鍵。 6. 欲清除儲存的基本重量，常按 <b>Mode</b> 至顯示 Count 模式後螢幕顯示 CLR.RPU 時按 <b>Yes</b> 確認</p>	 <p>The screenshot shows the NAVIGATOR scale display with '1300 g' on the screen. Below the screen are four buttons: 'Zero' (with a circled '1' above it), 'Print Units', 'Function Mode', and 'Tare Menu'. Under 'Zero' are 'Yes' and 'No' options. Under 'Print Units' is 'No'. Under 'Function Mode' is 'Back'. Under 'Tare Menu' is 'Exit'. A hand icon is pointing to the 'Function Mode' button.</p>

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